



## **OptiMICE Bottle Cap Tool (P/N C79190)**

Quick, ergonomic removal of bottle caps can be done with the bottle cap tool included with every rack that uses water bottles. The tool shown in (Fig. 1) is shipped, ready for installation on a counter or table. The tool can also be used un-secured after installing the flexible handle grip (Fig. 2).



Fig. 1



Fig. 2

### **Using the tool un-secured:**

1. Install the flexible handle grip.
2. Place the tool under the bottle cap with the tabs pointing down as shown in Fig. 3. Place index finger on the bottle cap to secure (Fig. 4) and lift the cap off.

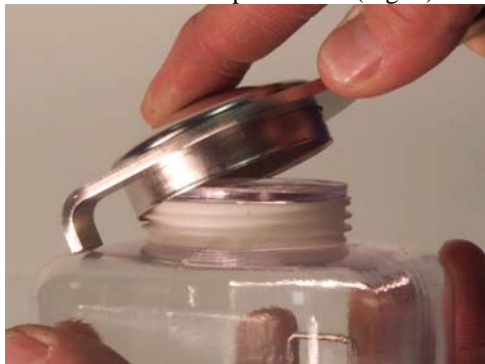


Fig. 3



Fig. 4

### **Using the tool secured to a counter or table:**

1. Mount the tool to the counter using the two holes pre-drilled for you.
2. Place the bottle cap in the tool and pull the bottle down and towards you (Fig. 5).



Fig. 5