



**ANIMAL CARE
SYSTEMS**

7086 South Revere Parkway, Suite #100, Centennial, Colorado 80112

Toll Free 888-827-3861 Phone 720-283-0177 Fax 720-283-0179

www.**AnimalCareSystems.com**

Cirrus SPF® Misting Station UNPACKING INSTRUCTIONS:

USE CAUTION WHEN LIFTING!

Pallet is accessible from two ends with a forklift or pallet jack. Be sure your equipment can handle the weight and size. Weights and dimensions are noted below.

TOOLS REQUIRED: Phillips driver and cutters or utility knife.

1. Remove banding with box cutters.
2. Remove the tape on the top of the box with box cutters.
3. Remove the cardboard panel.
4. Remove the banding holding the unit to the wooden pallet.
5. With 2 people, carefully rock the misting station up onto its casters, keeping the base against the pallet.



Remove banding with box cutters



Remove the tape with box cutters



Remove the cardboard panel.



Remove the banding holding the unit to the wooden pallet



Lift and rotate on to casters.



| In (mm) | Unit with Pallet & Box |
|---------------|------------------------|
| Length | 56" (142.25 cm) |
| Width | 32" (81.25 cm) |
| Height | 32" (81.25 cm) |
| Weight | 220 lbs. (99.79 kg) |

Copies of these instructions, and other useful information, are also available for download at:
www.animalcaresystems.com