



### OptiMICE® UNPACKING INSTRUCTIONS:

Copies of these instructions are also available for download in a PDF format at: [www.animalcaresystems.com](http://www.animalcaresystems.com)

Pallet is accessible from two sides with a forklift or pallet jack. Be sure your equipment can handle the weight and size. USE CAUTION WHEN LIFTING RACKS!! RACKS ARE TOP HEAVY!! Rack weights and dimensions are noted below.

**IF YOUR RACKS ARE SHIPPED VIA OCEAN CONTAINER (NOT BOXED) PLEASE SKIP TO STEP #5**

**TOOLS REQUIRED: Phillips driver, strap cutters and 7/16" (11mm) wrench.**

1. Cut the vertical and horizontal straps.
2. Remove the cardboard top.
3. Remove the two large cardboard sides.
4. Remove the 4 cardboard/foam corners from the rack.

**CAUTION: Check the corner pieces for any accessories that may be packed inside! They will be marked with yellow CAUTION TAPE.**




5. Cut the two plastic tie wraps holding the casters in place.




OptiMICE® Racks		
	Rack Only	Rack/pallet & box
	Lbs (Kg)	Lbs (Kg)
Standard Rack	560 (254)	590 (268)
Rack w/auto water	586 (266)	616 (279)
	Inches (mm)	Inches (mm)
Length	40 (1016)	44 (1118)
Width	34 (864)	44 (1118)
Height	78 (1981)	79 (2006)

6. Remove all four UPPER hex bolts (connecting the L-brackets to the rack) with 7/16" (or 11mm) wrench.

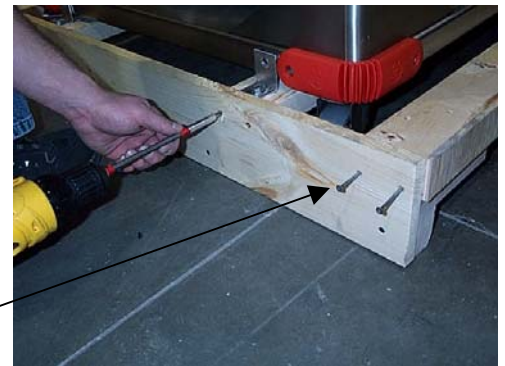

7. Remove the eight (8) screws from the 1x6 on one side using a Phillips driver.

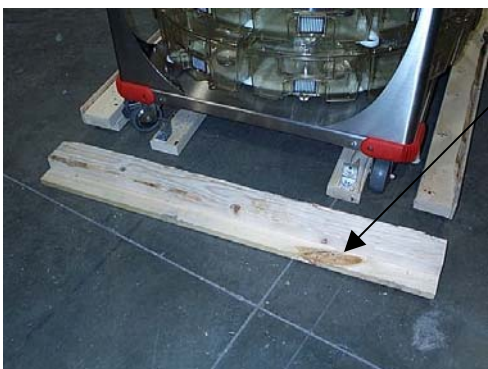

8. Knock off the 1x6 end piece. Unlock the two casters.



10. Roll the rack off the remaining pallet.