



OptiMICE C79072 CAGE FILTER REPLACEMENT

Old Filter Removal:

From the outside of the cage push the filter out with both thumbs (Figure 1). Repeat on the opposite end of the cage.



Figure 1: Removing Filter

Filter Installation:

1. Moisten the seals on the filter assemblies with weak soapy water.
2. From inside the cage, place the filter in the opening so that the rib on the seal is against the cage all around (Figures 2 and 3).

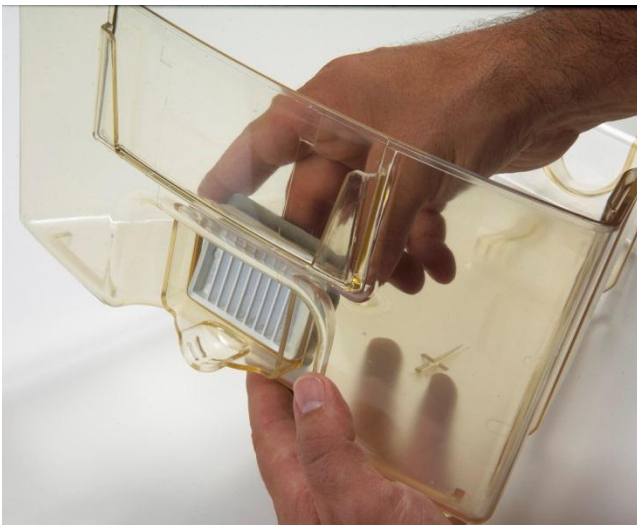


Figure 2: Installing Filter



Figure 3: Installing Filter



3. Push the top portion of the filter into the cage until it snaps into place (Figure 4). Push the bottom portion of the filter into the cage until it snaps into place (Figure 5).

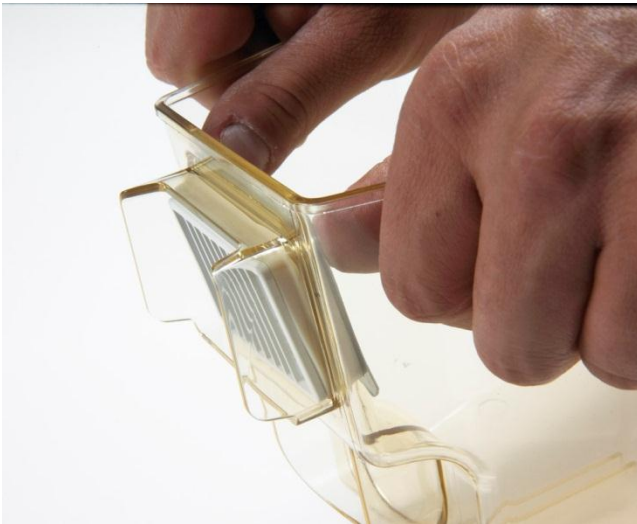


Figure 4: Installing Top Portion

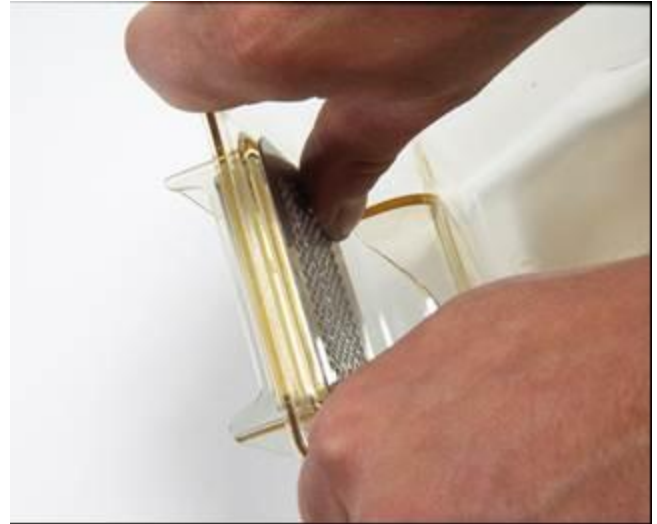


Figure 5: Installing Bottom Portion

(NOTE: it is important to push the filter into the cage in a two-step process to avoid damage to the seal.)

Filter Inspection:

4. Inspect the seal and be sure the ribs are straight and the seal has no tears (Figure 6).

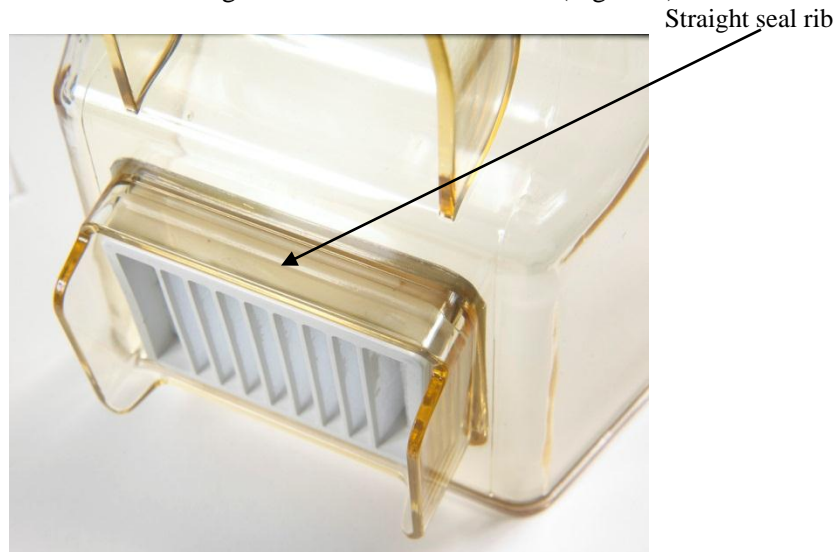


Figure 6: Inspecting the Seal