



ANIMAL CARE  
SYSTEMS

7086 South Revere Parkway, Suite #100, Centennial, Colorado 80112

Toll Free 888-827-3861 Phone 720-283-0177 Fax 720-283-0179

www.**AnimalCareSystems.com**

### OptiMICE® UNPACKING INSTRUCTIONS:

**USE CAUTION WHEN LIFTING! RACKS ARE TOP HEAVY!**

Pallet is accessible from two sides with a forklift or pallet jack. Be sure your equipment can handle the weight and size. Rack weights and dimensions are noted below.

**TOOLS REQUIRED: Phillips driver, cutters or utility knife, and 7/16" (11mm) wrench.**

(Racks shipped via ocean container will not be boxed. Please skip to step 4.)

1. Cut the vertical and horizontal straps.
2. Remove the cardboard top.
3. Remove the two large cardboard sides.
4. Remove the 4 cardboard/foam corners from the rack.
5. Carefully cut and remove the protective film wrapping.

*Check the corner pieces for any accessories that may be packed inside. They will be labeled and/or marked with yellow caution tape.*



Cut the two plastic ties attached to the front locking casters.



Remove all four (4) L-brackets holding rack to pallet with a 7/16" (or 11mm) wrench.



Remove ten (10) wood screws from the top of the pallet using a Phillips screwdriver.



Remove screws (two in front and back) from center support and remove 2 x 4.



Knock off the 1 x 6 end piece of the pallet and unlock the two front casters.



Roll the rack off of the remaining pallet.

Lbs (Kg)	Rack Only	Rack with Pallet & Box
<b>Standard Rack</b>	560 (254)	590 (268)
<b>Rack with Auto Water</b>	586 (266)	616 (279)

In (mm)	Rack Only	Rack with Pallet & Box
<b>Length</b>	40 (1016)	44 (1118)
<b>Width</b>	34 (864)	44 (1118)
<b>Height</b>	78 (1981)	79 (2006)

Copies of these instructions, and other useful information, are also available for download at: [www.animalcaresystems.com](http://www.animalcaresystems.com)